

ApoAlert™ DNA Fragmentation Assay Kit User Manual



Clontech

United States/Canada
800.662.2566

Asia Pacific
+1.650.919.7300

Europe
+33.(0)1.3904.6880

Japan
+81.(0)77.543.6116

Clontech Laboratories, Inc.
A Takara Bio Company
1290 Terra Bella Ave.
Mountain View, CA 94043
Technical Support (US)
E-mail: tech@clontech.com
www.clontech.com

Cat. Nos. 630107
630108
PT3137-1 (PR143851)
Published April 2011

Table of Contents

I. Introduction	3
II. List of Components	5
III. Additional Materials Required	6
IV. Controls	8
V. Sample Preparation for Microscopic Detection	9
A. Adherent Cells	9
B. Suspended Cell Cultures	10
C. Tissue Sections	11
VI. Detection by Microscopy	13
VII. Detection by Flow Cytometry	15
VIII. Troubleshooting Guide	17
IX. References	18

List of Tables

Table I: Modes of Action of Apoptosis-Inducing Reagents	8
---	---

Contact Us For Assistance

Customer Service/Ordering:	Technical Support:
Telephone: 800.662.2566 (toll-free)	Telephone: 800.662.2566 (toll-free)
Fax: 800.424.1350 (toll-free)	Fax: 650.424.1064
Web: www.clontech.com	Web: www.clontech.com
E-mail: orders@clontech.com	E-mail: tech@clontech.com

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product's webpage at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

Clontech, the Clontech logo, ApoAlert, and Living Colors are trademarks of Clontech Laboratories, Inc. All other marks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech is a Takara Bio Company. ©2011 Clontech Laboratories, Inc.

I. Introduction

Apoptosis, or programmed cell death, plays a fundamental role in many normal biological processes, as well as in several disease states (Wyllie *et al.*, 1980; Ellis *et al.*, 1991; Cohen *et al.*, 1992; Thompson, 1995; Nicholson, 1996). Apoptotic cells undergo many distinct morphological and biochemical changes, including fragmentation of nuclear DNA (Schwartzman & Cidlowski, 1993; Walker *et al.*, 1991; Oberhammer *et al.*, 1993; Roy *et al.*, 1992). During apoptosis, cellular endonucleases cleave nuclear DNA between nucleosomes, producing a mixture of DNA fragments whose lengths vary in multiples of 180 to 200 bp.

The ApoAlert DNA Fragmentation Assay Kit detects apoptosis-induced nuclear DNA fragmentation via a fluorescence assay. The assay is based on terminal deoxynucleotidyl transferase (TdT)-mediated dUTP nick-end-labeling (TUNEL; Gavrieli *et al.*, 1992; Piqueras *et al.*, 1996; Facchinetti *et al.*, 1991; Wijsman *et al.*, 1993). TdT catalyzes incorporation of fluorescein-dUTP at the free 3'-hydroxyl ends of fragmented DNA; fluorescein-labeled DNA can be detected via fluorescence microscopy or flow cytometry. This kit uses a direct labeling procedure for detecting DNA fragmentation at the single-cell level. With a fluorescence microscope equipped with FITC filters, you can visualize areas of apoptotic cells in situ in tissue sections. The kit can be used with adherent cells and cells in suspension.

Other ApoAlert Apoptosis Detection Kits

Although the TUNEL assay is widely used for detecting apoptosis, DNA fragmentation occurs relatively late in the process. Recent studies have elucidated some molecular events that are characteristic of the early stages of apoptosis (Zhang, *et al.*, 1996; Koester, *et al.*, 1997; Green, 1998; Nicholson & Thornberry, 1997; Thornberry & Littlewood, 1998). As a result, the TUNEL assay can be complemented with other assays that detect apoptosis at earlier stages.

- The **ApoAlert Caspase-3 Fluorescent Assay Kits** (Cat. Nos. 630214 & 630215) and **ApoAlert Caspase-3 Colorimetric Assay Kits** (Cat. Nos. 630216 & 630217) provide assays for caspase-3 protease activity, or other proteases that recognize the caspase-3 cleavage sequence, DEVD. Induction of caspase-3 activity is one of the earliest events in Fas receptor-induced apoptosis; however, caspase-3 activity is induced somewhat later in staurosporine-induced apoptosis.
- The **ApoAlert Caspase-8 Fluorescent Assay Kits** (Cat. Nos. 630218 & 630219) and **ApoAlert Caspase-8 Colorimetric Assay Kits** (Cat. Nos. 630220 & 630221) provide assays for caspase-8 protease activity. Induction of caspase-8 is one of the earliest events in Fas receptor-induced apoptosis. Note that staurosporine does not induce caspase-8.

I. Introduction *continued*

- The **ApoAlert Annexin V Apoptosis Kit (FITC)**: [Cat. Nos. 630109 & 630110] is based on externalization of phosphatidylserine (PS) from the inner to the outer leaflet of the plasma membrane. This event occurs shortly after caspase activation. Annexin V is a protein with strong affinity for PS. The assay is simple and can be used with fluorescence microscopy or flow cytometry.
- The **ApoAlert Cell Fractionation Kit** (Cat. No. 630105) provides a method for isolating a mitochondria-enriched fraction from the cytoplasm of apoptotic and non-apoptotic cells. It includes two different antibodies against Cytochrome C. The first antibody is used to identify the mitochondria-enriched fraction. The second antibody is used to determine if Cytochrome C has been released from the mitochondria and is present in the cytosolic fraction—an indicator of mitochondrial involvement in apoptosis.
- The **ApoAlert Caspase-3 Assay Plate** (Cat. No. 630223) is a 96-well plate with the fluorogenic substrate of caspase-3 immobilized in the wells. When cell lysate containing active caspase-3 is placed in the wells, caspase-3 cleaves the substrate, releasing a fluorescent product that can be detected with a standard absorbance microplate reader.
- The **ApoAlert Caspase Profiling Plate** (Cat. Nos. 630225 & 630226) is a 96-well plate with fluorogenic substrates for four different caspases immobilized in separate wells. When cell lysate containing the active caspase of interest is added to the wells, the caspase cleaves the substrate, yielding a fluorescent product that can be detected with a standard fluorescence microplate reader.
- The **ApoAlert pDsRed2-Bid Vector** (Cat. No. 632419) encodes a fusion of a Living Colors® Fluorescent Protein (DsRed2), a red fluorescent protein, and Bid, a member of the Bcl-2 family. In healthy, non-apoptotic cells, Bid resides in the cytosol as soluble protein. Upon induction of apoptosis, Bid trafficks to the mitochondria. In cells expressing the Bid-DsRed fusion, Bid trafficking can be detected by fluorescence microscopy.

Clontech also offers apoptosis-inducing reagents as well as antibodies, protease inhibitors, stains, and other reagents for apoptosis research.

II. List of Components

Upon receipt, store the kit at -20°C .

After first use, store 20X SSC and plastic coverslips at room temperature. Store all other components at -20°C .

Important: The nucleotide mix is sensitive to light.

	Cat. No. 630107 (25 assays)	Cat. No. 630108 (100 assays)
• Equilibration Buffer	4 x 1 ml	10 x 1.6 ml
• Nucleotide Mix	125 μl	500 μl
• TdT Terminal Transferase (45 U/ μl)	28 μl	100 μl
• 20X SSC (see note below)	29 ml	2 x 58 ml
• Proteinase K (10 mg/ml)	440 μl	1.85 ml
• Plastic Coverslips	25	100

Composition of components

- Equilibration Buffer:

200 mM	Potassium Cacodylate (pH 6.6 at 25°C)
25 mM	Tris HCl (pH 6.6 at 25°C)
0.2 mM	DTT
0.25 mg/ml	BSA
2.5 mM	Cobalt chloride

Potassium Cacodylate (dimethylarsinic acid) is a mild irritant. Avoid skin and eye contact by wearing gloves and safety glasses.

- 20X SSC:

3 M	NaCl
300 mM	$\text{Na}_3\text{Citrate} \bullet \text{H}_2\text{O}$

If stored at -20°C , 20X SSC must be warmed to room temperature to ensure that all salts are in solution prior to 10X dilution with deionized H_2O . After opening the kit, 20X SSC can be stored at room temperature.

III. Additional Materials Required

For working with cell cultures (Sections V.A & V.B)

- Poly-L-lysine coated microscope slides

Apply 50–100 µl of a 0.1% (w/v) aqueous solution of poly-L-lysine (Sigma Cat. No. P8920) to the surface of sterile glass microscope slide. Spread evenly over the portion of the slide that will be used for growing and fixing cells. Air-dry the slides for 1 hr and store at 4°C for up to one week. Alternatively, use Lab-Tek Chamber Slides or equivalent commercial slides.

- Coplin jars filled with:

- Phosphate-buffered saline (PBS [pH 7.4]; 6 jars)
- 4% Formaldehyde/PBS (pH 7.4; 1 jar)

Prepare formaldehyde/PBS solutions immediately before use. Combine 70 ml of PBS and 25 ml of 16% methanol-free formaldehyde (Polyscience, Inc. Cat. No. 18814 or Aldrich Cat. No. 44124-4). Adjust the pH to 7.4 with 1 N NaOH.

- 0.2% Triton X-100/PBS (1 jar)
- Deionized H₂O (1 jar; reserve for DNase-treated samples)
- DNase I Buffer

40 mM	Tris-HCl (pH 7.9)
10 mM	NaCl
6 mM	MgCl ₂
10 mM	CaCl ₂

- DNase I (0.5 µg to 1 µg/ml = final concentration)
- [Suspension cells only] Refrigerated centrifuge for collecting cells

For working with paraffin-embedded tissue sections (Section V.C)

- Coplin jars filled with:

- Xylene (2 jars)
- Ethanol rehydration series:
 - 100% ethanol (2 jars)
 - 95% ethanol (1 jar)
 - 85% ethanol (1 jar)
 - 70% ethanol (1 jar)
 - 50% ethanol (1 jar)

- 0.85% NaCl (1 jar)
- PBS (7 jars)
- 4% Formaldehyde/PBS (pH 7.4) (2 jars)
- 20 µg/ml Proteinase K solution (1 jar)

Combine 2 µl of 10 mg/ml Proteinase K (provided in the kit) and 998 µl of 100 mM Tris-HCl (pH 8.0), 50 mM EDTA.

III. Additional Materials Required *continued*

For microscopic detection of apoptosis (Section VI)

- Forceps for handling slides and coverslips
- 37°C incubator (dark, humidified)
- 2X SSC
Dilute 20X SSC (included in the kit) 10-fold with H₂O.
- Coplin jars filled with
 - PBS (3 jars)
 - Propidium iodide/PBS (PI/PBS; 1 jar), 10 µg/ml = Final Concentration.
Stock: Dissolve 10 mg of propidium iodide (Sigma, Cat. No. P4170) in 10 ml of PBS. Store in the dark at 4°C.
 - Deionized H₂O (3 jars)
- PI/RNase/PBS
 - 10 µg/ml Propidium iodide
 - 0.5 µg/ml DNase-free RNase
- Anti-Fade reagent (Molecular Probes, Cat. No. S7461)

For detection of apoptosis by flow cytometry (Section VII)

- Refrigerated centrifuge for collecting cells
- 37°C water bath
- PBS
- 1% Formaldehyde/PBS (pH 7.4)
Prepare formaldehyde/PBS solutions immediately before use. Combine 90 ml of PBS and 6.25 ml of 16% methanol-free formaldehyde. Adjust the pH to 7.4 with 1 N NaOH.
- 20 mM EDTA (pH 8.0)
- 0.1% Triton X-100/BSA/PBS
Combine:

5 ml	PBS
5 ml	0.2% Triton X-100
50 mg	BSA
- Propidium iodide/PBS (PI/PBS; 1 jar)
Dissolve 10 mg of propidium iodide (Sigma, Cat. No. P4170) in 10 ml of PBS. Store in the dark at 4°C.
- PI/RNase/PBS
 - 2.5 µg/ml Propidium iodide
 - 0.5 µg/ml DNase-free RNase

IV. Controls

We recommend including positive and negative controls.

There are two types of positive controls:

- **Biological controls.** Use an apoptosis-inducing reagent to induce apoptosis. Table I describes the modes of action of several such reagents; some are available from Clontech.

TABLE I: MODES OF ACTION OF APOPTOSIS-INDUCING REAGENTS

<u>Agent</u>	<u>Mode of Action</u>
Actinomycin D	Inhibits RNA synthesis
Anti-Fas antibody (clone CH-11)	Binds and activates Fas cell-surface protein (CD95)
Cycloheximide	Inhibits protein synthesis
C2-Ceramide	Second messenger in sphingomyelin pathway
Dexamethasone	Binds glucocorticoid receptor
Etoposide	Inhibits topoisomerase
Human TNF- α	Binds TNF receptor when applied with cyclohexamide
Staurosporine	Inhibits protein kinases (broad spectrum)
Vincristine Sulfate	Disrupts microtubules

- **DNase-treated controls.** Prepare as described in Step V.A.8. DNase I treatment creates free 3'-OH ends that are labeled with fluorescein dUTP. Typically, >70% of treated cells will give a positive fluorescent signal.

There are two types of negative controls:

- **Biological controls.** Cells or tissues are not subjected to apoptosis-inducing treatments.
- **TdT-minus controls.** Cells or tissues are not treated with TdT.

V. Sample Preparation for Microscopic Detection

A. Adherent Cells

Most steps are performed in Coplin jars (Section III). **Keep Coplin jars that have been used with DNase separate from DNase-free Coplin jars.**

1. Grow cells on Lab-Tek Chamber Slides, equivalent commercial slides, or on poly-L-lysine-coated slides prepared in your laboratory (see Section III).
2. Induce apoptosis in samples.
3. Following induction, wash the slides by dipping twice in a Coplin jar containing PBS. Use a fresh jar for each wash.
4. To fix the cells, immerse the slides in a Coplin jar containing fresh 4% formaldehyde/PBS at 4°C for 25 min.
Note: You may store fixed cells for up to two weeks at -20°C in 70% ethanol.
5. Immerse the slides in a Coplin jar containing fresh PBS for 5 min at room temperature.
6. Repeat PBS wash.
7. Permeabilize the cells by immersing the slides in a Coplin jar containing prechilled 0.2% Triton X-100/PBS. Incubate for 5 min on ice.
8. [Optional] Prepare a DNase-treated positive control:
 - a. Add 100 µl of DNase I Buffer.
 - b. Incubate at room temperature for 5 min.
 - c. Gently tap the slide to remove the liquid.
 - d. Add 100 µl of DNase I Buffer containing 0.5–1 µg/ml DNase.
 - e. Incubate at room temperature for 10 min.
 - f. Gently tap the slide to remove the liquid.
 - g. Immerse 3–4 times in H₂O in a Coplin jar reserved for DNase-treated samples.
9. Immerse the slides in a Coplin jar containing fresh PBS for 5 min at room temperature.
10. Transfer the slides to a second Coplin jar containing fresh PBS for 5 min at room temperature.
11. Proceed to Section VI.

V. Sample Preparation for Microscopic Detection *continued*

B. Cultures of Suspended Cells

1. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
2. Aspirate the supernatant carefully.
3. Add 500 µl of PBS and gently resuspend the cells.
4. Centrifuge the cells at 300 x g for 10 min at 4°C.
5. Aspirate the supernatant carefully.
6. Resuspend in PBS at a concentration of 2×10^7 cells/ml.
7. Transfer 50–100 µl of the cell suspension onto a poly-L-lysine-coated or silanized slide.
8. GENTLY spread the cell suspension with a clean slide.
9. Fix cells by immersing the slides in a Coplin jar containing fresh 4% formaldehyde/PBS at 4°C for 25 min.
Note: You may store fixed cells for up to two weeks at –20°C in 70% ethanol.
10. Immerse the slides in a Coplin jar containing fresh PBS for 5 min at room temperature.
11. Repeat PBS wash.
12. Permeabilize the cells by immersing the slides in a Coplin jar containing prechilled 0.2% Triton X-100/PBS. Incubate 5 min on ice.
13. [Optional] Prepare a DNase-treated positive control:
 - a. Add 100 µl of DNase I Buffer.
 - b. Incubate at room temperature for 5 min.
 - c. Gently tap the slide to remove the liquid.
 - d. Add 100 µl of DNase I Buffer containing 0.5–1 µg/ml DNase.
 - e. Incubate at room temperature for 10 min.
 - f. Gently tap the slide to remove the liquid.
 - g. Immerse 3–4 times in H₂O in a Coplin jar reserved for DNase-treated samples.
14. Immerse the slides in a Coplin jar containing fresh PBS for 5 min at room temperature.
15. Transfer the slides to a second Coplin jar containing fresh PBS for 5 min at room temperature.
16. Proceed to Section VI.

V. Sample Preparation for Microscopic Detection *continued*

C. Tissue Sections

This protocol describes the preparation of formalin-fixed, paraffin-embedded tissue sections mounted on glass slides. For information on fixing and embedding techniques, see Ben-Sasson et al. (1995). Most steps are performed in Coplin jars (Section III). **You must separate Coplin jars that have been used with DNase from DNase-free Coplin jars.**

Note: If using fresh frozen tissue sections, proceed directly to Step 7.

1. Remove paraffin by immersing the slides in a Coplin jar containing fresh xylene. Incubate at room temperature for 5 min.
2. Repeat in a second Coplin jar containing fresh xylene.
3. Immerse the slides in a Coplin jar containing 100% ethanol and incubate at room temperature for 5 min.
4. Rehydrate the slides by sequential 3 min, room temperature incubations in Coplin jars containing:
 - 100% ethanol
 - 95% ethanol
 - 85% ethanol
 - 70% ethanol
 - 50% ethanol
5. Immerse the slides in a Coplin jar containing 0.85% NaCl and incubate at room temperature for 5 min.
6. Immerse the slides in a Coplin jar containing PBS and incubate at room temperature for 5 min.
7. Fix the slides by immersing them in a Coplin jar containing fresh 4% formaldehyde/PBS, and incubate at room temperature for 15 min.
8. Wash the slides by immersing them in a Coplin jar containing PBS, and incubate at room temperature for 5 min.
9. Transfer to another Coplin jar containing PBS, and incubate at room temperature for 5 min.
10. Allow the liquid to drain thoroughly and place the slides on a flat surface.
11. Prepare 20 µg/ml Proteinase K solution (Section III) and cover each section with 100 µl of it. Incubate at room temperature for 5 min.

Note: Optimizing the length of Proteinase K treatment may be useful. For example, sections thicker than 5 µm may require longer treatments. However, overtreatment may loosen sections, causing them to wash off the slides during subsequent steps.

V. Sample Preparation for Microscopic Detection *continued*

12. Immerse the slides in a Coplin jar containing PBS and incubate at room temperature for 5 min.
13. Transfer the slides to a Coplin jar containing 4% formaldehyde/PBS, and incubate at room temperature for 5 min.
14. Wash the slides by immersion in a Coplin jar containing PBS and incubate at room temperature for 5 min.
15. Proceed to Section VI.

VI. Detection by Microscopy

This assay procedure can be used for any sample obtained in Section V. See Section III for information about Coplin jars and solutions. **Separate Coplin jars that have been used with DNase from DNase-free Coplin jars.**

1. Remove the slides from PBS and tap gently to remove excess liquid.
2. Cover the cells in 100 μ l of equilibration buffer.
3. Using forceps, gently place a piece of plastic coverslip on top of the cells to evenly spread the liquid.
4. Equilibrate at room temperature for 10 min.
5. Thaw Nucleotide Mix on ice and prepare TdT incubation buffer for all experimental samples, positive controls, and biological negative controls.

Component	Volume	Total No. of Reactions	Total Volume
Equilibration Buffer	45 μ l	—	— μ l
Nucleotide Mix	5 μ l	—	— μ l
TdT Enzyme	1 μ l	—	— μ l
	51 μ l	—	— μ l

6. For your TdT-minus negative control, prepare TdT-minus control incubation buffer by replacing the TdT Enzyme with deionized H₂O.

Note: Protect Nucleotide Mix and TdT incubation buffer from light. Keep on ice at all times.

7. Using forceps, remove the plastic coverslip and gently tap the slides to remove excess liquid.
 8. Carefully blot dry around the edges with tissue paper.
 9. Gently place 50 μ l of TdT incubation buffer on the cells on a ~5 cm² area.
- Note:** Do not allow the cells to dry out.
10. [Optional] If performing a TdT-minus negative control, add 50 μ l of TdT-minus control incubation buffer.

From this point, protect samples from light by wrapping Coplin jars in aluminum foil or keeping the jars in a covered box.

11. Using forceps, gently place a piece of plastic coverslip on top of the cells to evenly spread the liquid.
12. To perform the tailing reaction, place the slides in a dark, humidified 37°C incubator for 60 min.

Note: Ensure high humidity by placing wet paper towels in the bottom of dry incubators.

VI. Detection by Microscopy *continued*

13. Using forceps, remove the plastic coverslips.
14. Terminate the tailing reaction by immersing the slides in a Coplin jar containing 2X SSC. Incubate at room temperature for 15 min.
15. Wash the cells by transferring the slides to a fresh Coplin jar filled with PBS. Incubate at room temperature for 5 min.
16. Repeat Step 15 twice.
17. Stain the cells with propidium iodide (PI) by incubating the slides in a Coplin jar containing 40 ml of fresh PI/PBS at room temperature for 5–10 min.
Note: If PI staining is too intense, include 0.5 µg/ml of DNase-free RNase at this step.
18. Wash the cells by transferring the slides to a fresh Coplin jar filled with deionized H₂O and incubate at room temperature for 5 min.
19. Repeat Step 18 twice.
20. [Optional] Add a drop of Anti-Fade reagent and cover the treated portion of the slide with a glass coverslip.
21. [Optional] Seal the edges of the coverslip with rubber cement or clear nail polish and allow to dry for at least 5–10 min.

You may store sealed slides overnight at 4°C in the dark.

View the slides as soon as possible. Apoptotic cells will exhibit strong, nuclear green fluorescence using a standard fluorescein filter set (520 ± 20 nm). All cells stained with PI exhibit strong red cytoplasmic fluorescence when viewed at >620 nm.

VII. Detection by Flow Cytometry

1. Collect $3-5 \times 10^6$ cells by centrifugation at 300 x g for 10 min at 4°C.
2. Aspirate the supernatant carefully.
3. Add 5 ml of PBS and gently resuspend the cells.
4. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
5. Aspirate the supernatant carefully.
6. Repeat Steps 3–5.
7. Resuspend the cells in 0.5 ml of PBS.
8. Fix the cells by adding 5 ml of fresh, prechilled 1% formaldehyde/PBS. Incubate at 4°C for 20 min.
9. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
10. Carefully aspirate the supernatant.
11. Gently resuspend the cells in 5 ml of PBS.
12. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
13. Carefully aspirate the supernatant.
14. Repeat steps 11–13.
15. Gently resuspend the cells in 0.5 ml of PBS.
16. Permeabilize the cells by adding 5 ml of 70% ice-cold ethanol. Incubate at –20°C for at least 4 hr. The cells can be stored at –20°C for up to 1 week.
Note: For a faster procedure, permeabilize the cells by incubating on ice with 5 ml of prechilled 0.2% Triton X-100.
17. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
18. Remove supernatant and gently resuspend the cells in 5 ml of PBS.
19. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
20. Remove the supernatant and gently resuspend the cells in 1 ml of PBS.
21. Transfer 2×10^6 cells to an amber 1.5-ml microcentrifuge tube.
Note: Protect samples from light from this point.
22. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
23. Carefully aspirate the PBS.
24. Gently resuspend the cells in 80 µl of Equilibration Buffer.
25. Incubate at room temperature for 5 min.

VII. Detection by Flow Cytometry *continued*

26. Thaw the Nucleotide Mix on ice and prepare TdT incubation buffer for all experimental samples, positive controls, and biological negative controls.

Component	Volume	Total No. of Reactions	Total Volume
Equilibration Buffer	45 μ l	—	— μ l
Nucleotide Mix	5 μ l	—	— μ l
TdT Enzyme	1 μ l	—	— μ l
	51 μ l	—	— μ l

27. For your TdT-minus negative control, prepare TdT-minus control incubation buffer by replacing the TdT Enzyme with deionized H₂O.

Note: Protect Nucleotide Mix and TdT incubation buffer from light. Keep on ice at all times.

28. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
29. Carefully aspirate the supernatant.
30. Resuspend in 50 μ l of TdT incubation buffer.
31. Incubate at 37°C in a water bath for 60 min, protected from direct light. Mix the cell suspension by gently pipetting every 15 min.
32. Add 1 ml of 20 mM EDTA to terminate the reaction, and mix by gentle vortexing.
33. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
34. Carefully aspirate the supernatant.
35. Resuspend the cells in 1 ml of 0.1% Triton X-100/BSA/PBS.
36. Repeat Steps 33–35.
37. Collect the cells by centrifugation at 300 x g for 10 min at 4°C.
38. Carefully aspirate the supernatant. Stain the cells with propidium iodide (PI) by gently resuspending them in 0.5 ml of PI/RNase/PBS.
39. Incubate at room temperature in the dark for 15–30 min.

Analyze the cells as soon as possible. Apoptotic cells can be collected based on green fluorescence at 520 $\pm$ 20 nm. All cells can be sorted based on red fluorescence at >620 nm.

VIII. Analysis of Results & Troubleshooting Guide

- **High background**

High background can be caused by nonspecific incorporation of fluorescein-12-dUTP. This problem can occur if the cells dry out at Step VI.11 or if they are insufficiently washed at Step VI.17. Correct for insufficient washing by washing the slides 3 x 5 min with Triton X-100/BSA/PBS.

- **No positive signal**

This problem usually indicates poor permeabilization with Triton X-100 (Steps V.A.7, V.B.12, or VII.16), or with Proteinase K (Step V.C.11). Address the problem by incubating longer with the permeabilization reagent.

- **Tissue sections are loose**

Tissue slides can loosen for two reasons: 1) poor preparation of the slides prior to mounting of tissue sections; or 2) overdigestion by Proteinase K (Step V.C.11). Correct the first problem by switching from poly-L-lysine-coated slides to slides coated with 3-aminopropyl triethoxysilane (Ben-Sasson et al. 1995). Correct the second problem by optimizing the Proteinase K incubation time.

- **Too few cells for analysis**

You can lose cells at several points. If this problem occurs consistently, first increase the starting number of cells. If you are preparing suspension cells for attachment to slides (Section V.B), try washing the cells with PBS + 1% BSA during the centrifugation steps. If a cytospin centrifuge is available, use it to prepare your samples. Finally, try washing the cells with PBS + 1% BSA during the centrifugation steps in Section VII.

IX. References

- Ben-Sasson, S. A., Sherman, Y. & Gavrieli, Y. (1995) Identification of dying cells by in situ staining. *Methods Cell. Biol.* **46**:29–39.
- Cohen, J. J., Duke, R. C., Fadok, V. A. & Sellins, K. S. (1992) Apoptosis and cell death in immunity. *Ann. Rev. Immunol.* **10**:267–293.
- Ellis, R. E., Yuan, J. & Horvitz, H. R. (1991) Mechanisms and function of cell death. *Ann. Rev. Cell Biol.* **7**:663–698.
- Facchinetti, A., Tessarollo, L., Mazzocchi, M., Kingston, R., Collavo, D. & Biasi, G. (1991) An improved method for the detection of DNA fragmentation. *J. Immunol. Methods* **136**:1251–1256.
- Gavrieli, Y., Sherman, Y. & Ben-Sasson, S. A. (1992) Identification of programmed cell death in situ via specific labeling of nuclear DNA fragmentation. *J. Cell. Biol.* **119**:493–501.
- Green, D. R. (1998) Apoptotic pathways: the roads to ruin. *Cell* **94**:695–698.
- Koester, S. K., Roth, P., Mikulka, W. R., Schlossman, S. F., Zhang, C. & Bolton, W. E. (1997) Monitoring early cellular responses in apoptosis is aided by the mitochondrial membrane protein-specific monoclonal antibody APO2.7. *Cytometry* **29**:306–312.
- Nicholson, D. W. (1996) ICE/CED3-like proteases as therapeutic targets for the control of inappropriate apoptosis. *Nature Biotechnol.* **14**:297–301.
- Nicholson, D. W. & Thornberry, N. A. (1997) Caspases: killer proteases. *Trends Biochem. Sci.* **22**:299–306.
- Oberhammer, F., Wilson, J. W., Dive, C., Morris, I. D., Hickman, J. A., Wakeling, A. E., Walker, P. R., & Sikorska, M. (1993) Apoptotic death in epithelial cells: cleavage of DNA to 300 and/or 50 kb fragments prior to or in the absence of internucleosomal fragmentation. *EMBO J.* **12**:3679–3684.
- Piqueras, B., Autran, B., Debre, P., & Gorochoff, G. (1996) Detection of apoptosis at the single-cell level by direct incorporation of fluorescein-dUTP in DNA strand breaks. *BioTechniques* **20**:634–640.
- Roy, C., Brown, D. L., Little, J. E., Valentine, B. K., Walker, P. R., Sikorska, M., Leblanc, J., & Chaly, N. (1992) The topoisomerase II inhibitor teniposide (VM-26) induces apoptosis in unstimulated mature murine lymphocytes. *Exp. Cell. Res.* **200**:416–424.
- Schwartzman, R. A. & Cidlowski, J. A. (1993) Apoptosis: the biochemistry and molecular biology of programmed cell death. *Endocrine Rev.* **14**:133–151.
- Thompson, C. B. (1995) Apoptosis in the pathogenesis and treatment of disease. *Science* **267**:1456–1462.
- Thornberry, N. A. & Littlewood, Y. (1998) Caspases: Enemies Within. *Science* **281**:1312–1316.
- Walker, P. R., Smith, C., Youdale, T., Leblanc, J., Whitfield, J. F. & Sikorska, M. (1991) Topoisomerase II-reactive chemotherapeutic drugs induce apoptosis in thymocytes. *Cancer Res.* **51**: 1078–1085.
- Wijmsman, J. H., Jonker, R. R., Keijzer, R., Van de Velde, C. J. H., Cornelisse, C. J. & Van Dierendonck, J. H. (1993) A new method to detect apoptosis in paraffin sections: In situ end-labeling of fragmented DNA. *J. Histochem. Cytochem.* **41**:7–12.
- Wyllie, A. H., Kerr, J. F. R. & Currie, A. R. (1980) Cell death: The significance of apoptosis. *Intl. Rev. Cytol.* **68**:251–306.
- Zhang, C., Ao, Z., Seth A., & Schlossman, S. F. (1996) A mitochondrial membrane protein defined by a novel monoclonal antibody is preferentially detected in apoptotic cells. *J. Immunol.* **157**:3980–3987.