

**PRODUCT: pCMV-Myc & pCMV-HA Vector Set****CATALOG No.** 631604**LOT NUMBER**

Specified on product label.

**STORAGE CONDITIONS**

Store at -20°C

**SHELF LIFE**

1 year from date of receipt under proper storage conditions.

**SHIPPING CONDITIONS**

Dry ice (-70°C)

**DESCRIPTION**

A kit for expression, detection and coimmunoprecipitation of epitope-tagged proteins in mammalian cells. These vectors allow you to easily clone and express a gene of interest fused to an epitope tag. The antibodies to the tags allow detection and purification of the expressed tagged proteins.

**PACKAGE CONTENTS**

- 10 µg pCMV-Myc-N Vector (500 ng/µl)
- 10 µg pCMV-HA-N Vector (500 ng/µl)
- 500 µl c-Myc Monoclonal Antibody (200 µg/ml)
- 100 µl HA-Tag Polyclonal Antibody (IgG; 1 mg/ml)
- 50 µl pCMV Sequencing Primer (20 µM)

**SUGGESTED ANTIBODY DILUTIONS FOR WESTERN BLOTS:**

- c-Myc Monoclonal Antibody (1:100)
- HA-Tag Polyclonal Antibody (1:500)

**PRIMER SEQUENCE**

5'-GAT-CCG-GTA-CTA-GAG-GAA-CTG-AAA-AAC-3'  
(primer location indicated on Vector Information Packets)

**ANTIBODY SOURCES**

- c-Myc Monoclonal Antibody was raised in mice against a synthetic peptide (residues 408–439 of the human p62-c-Myc protein).
- HA-Tag Polyclonal Antibody was raised in rabbits against KLH-conjugated synthetic HA epitope (YPYDVPDYA).

**PRODUCT USER MANUALS**

User manuals for Clontech products are available for download at [www.clontech.com/manuals](http://www.clontech.com/manuals). The following user manuals apply to this product:

- pCMV-Myc-N Vector Information
- pCMV-HA-N Vector Information

**Clontech**

United States/Canada  
800.662.2566

Asia Pacific  
+1.650.919.7300

Europe  
+33.(0)1.3904.6880

Japan  
+81.(0)77.543.6116

Clontech Laboratories, Inc.  
A Takara Bio Company  
1290 Terra Bella Ave.  
Mountain View, CA 94043  
Technical Support (US)  
E-mail: [tech@clontech.com](mailto:tech@clontech.com)  
[www.clontech.com](http://www.clontech.com)

**FOR RESEARCH USE ONLY****QUALITY CONTROL DATA**

See back of page.

(110911)

**QUALITY CONTROL DATA****1. Plasmid identity and purity**

The identity of each plasmid was verified by: a) electrophoresis on an agarose/EtBr gel after digestion with the indicated enzymes, and b) sequence analysis of the epitope tag region. The purity of each plasmid was checked by determining the  $A_{260}/A_{280}$ .

<u>Vector</u>	<u>Enzyme</u>	<u>Fragment Size (kb)</u>
pCMV-Myc-N	EcoR I	3.8
	Pst I	2.7 & 1.1
pCMV-HA-N	EcoR I	3.8
	Pst I	2.7 & 1.1

**2. Antibodies**

The c-Myc and HA-Tag antibodies were tested by Western blotting. Cell lysates containing c-Myc- or HA-tagged proteins were electrophoresed on SDS polyacrylamide gels, followed by transfer to PVDF membranes. The blots were probed with 2 µg/ml c-Myc Monoclonal Antibody or 0.5 µg/ml HA-Tag Polyclonal Antibody, followed by secondary antibody.

## pCMV-Myc & pCMV-HA Vector Set

### CATALOG NO.

631604

### NOTICE TO PURCHASER:

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, in vitro diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without prior written approval of Clontech Laboratories, Inc.

Your use of this product is also subject to compliance with the licensing requirements listed below and described on the product's web page at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by these statements.

### TRADEMARKS:

Clontech and the Clontech logo are trademarks of Clontech Laboratories, Inc.

All other marks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech is a Takara Bio Company. ©2011 Clontech Laboratories, Inc. This document has been reviewed and approved by the Clontech Quality Assurance Department.

---

#### Clontech Laboratories, Inc.

A Takara Bio Company

1290 Terra Bella Avenue, Mountain View, CA 94043, USA

U.S. Technical Support: [tech@clontech.com](mailto:tech@clontech.com)

11/11/2011

#### United States/Canada

800.662.2566

#### Asia Pacific

+1.650.919.7300

#### Europe

+33.(0)1.3904.6880

#### Japan

+81.(0)77.543.6116