



pLVX-IRES-ZsGreen1 Vector Map and Multiple Cloning Site (MCS).

Description

pLVX-IRES-ZsGreen1 is an HIV-1-based, lentiviral expression vector that allows the simultaneous expression of your protein of interest and ZsGreen1 in virtually any mammalian cell type, including primary cells. ZsGreen1 is a human codon-optimized variant of the reef coral *Zoanthus sp.* green fluorescent protein, ZsGreen. The vector expresses the two proteins from a bicistronic mRNA transcript, allowing ZsGreen1 to be used as an indicator of transduction efficiency and a marker for selection by flow cytometry.

Expression of the bicistronic transcript is driven by the constitutively active human cytomegalovirus immediate early promoter ($P_{CMV IE}$) located just upstream of the MCS. An encephalomyocarditis virus (EMCV) internal ribosome entry site (IRES), positioned between the MCS and ZsGreen1, facilitates cap-independent translation of ZsGreen1 from an internal start site at the IRES/ZsGreen1 junction (1).

pLVX-IRES-ZsGreen1 contains all of the viral processing elements necessary for the production of replication-incompetent lentivirus, as well as elements to improve viral titer, transgene expression, and overall vector function. The woodchuck hepatitis virus posttranscriptional regulatory element (WPRE) promotes RNA processing events and enhances nuclear export of viral RNA (2), leading to increased viral titers from packaging cells. In addition, the vector includes a Rev-response element (RRE), which further increases viral titers by enhancing the transport of unspliced viral RNA out of the nucleus (3). Finally, pLVX-IRES-ZsGreen1 also contains a central polypurine tract/central termination sequence element (cPPT/CTS). During target cell infection, this element creates a central DNA flap that increases nuclear import of the viral genome, resulting in improved vector integration and more efficient transduction (4). The vector also contains a pUC origin of replication and an *E. coli* ampicillin resistance gene (Amp^r) for propagation and selection in bacteria.

(PR073579; published 13 July 2010)



Clontech

United States/Canada
800.662.2566

Asia Pacific
+1.650.919.7300

Europe
+33.(0)1.3904.6880

Japan
+81.(0)77.543.6116

Clontech Laboratories, Inc.
A Takara Bio Company
1290 Terra Bella Ave.
Mountain View, CA 94043
Technical Support (US)
E-mail: tech@clontech.com
www.clontech.com

Use

pLVX-IRES-ZsGreen1 is designed to constitutively coexpress your protein of interest and ZsGreen1 from P_{CMVIE} when transduced into mammalian cells. Before it can be transduced into target cells, the vector must be packaged into viral particles in HEK293T cells, using our Lenti-X™ HTX Packaging System (Cat. Nos. 631247 and 631249). This packaging system allows the safe production of high titer, infectious, replication-incompetent, VSV-G pseudotyped lentiviral particles that can infect a wide range of cell types, including nondividing and primary cells (5).

ZsGreen1 is the brightest commercially available green fluorescent protein. The presence of this protein allows transductants to be visualized by fluorescence microscopy and sorted by flow cytometry with standard FITC filter sets (ZsGreen1 has an excitation maximum of 493 nm and an emission maximum of 505 nm).

Location of Features

- 5' LTR (5' long terminal repeat): 1–635
- PBS (primer binding site): 636–653
- Ψ (packaging signal): 685–822
- RRE (Rev-response element): 1303–1536
- cPPT/CTS (central polypurine tract/central termination sequence): 2028–2151
- P_{CMVIE} (human cytomegalovirus immediate early promoter): 2185–2787
- MCS (multiple cloning site): 2803–2840
- IRES (encephalomyocarditis virus internal ribosome entry site): 2842–3416
- ZsGreen1 (*Zoanthus sp.* green fluorescent protein): 3417–4112
- WPRE (woodchuck hepatitis virus posttranscriptional regulatory element): 4126–4717
- 3' LTR (3' long terminal repeat): 4921–5557
- pUC origin of replication: 6027–6697 (complementary)
- Amp^r (ampicillin resistance gene; β -lactamase): 6842–7838 (complementary)

Selection of Transductants

- Marker: ZsGreen1

Propagation in *E. coli*

- Suitable host strains: DH5 α ™, DH10B™ and other general purpose strains.
- Selectable marker: plasmid confers resistance to ampicillin (100 μ g/ml) in *E. coli* hosts.
- *E. coli* replication origin: pUC
- Copy number: high

Notes:

The vector sequence was compiled from information in the sequence databases, published literature, and other sources, together with partial sequences obtained by Clontech. This vector has not been completely sequenced.

The viral supernatants produced by this lentiviral vector could contain potentially hazardous recombinant virus. Due caution must be exercised in the production and handling of recombinant lentivirus. Appropriate NIH, regional, and institutional guidelines apply.

References

1. Jang, S. K. *et al.* (1988) *J. Virol.* **62**(8):2636–2643.
2. Zufferey, R. *et al.* (1999) *J. Virol.* **73**(4):2886–2892.
3. Cochrane, A. W. *et al.* (1990) *Proc. Natl. Acad. Sci. USA* **87**(3):1198–1202.
4. Zennou, V. *et al.* (2000) *Cell* **101**(2):173–185.
5. Wu, X. *et al.* (2000) *Mol. Ther.* **2**(1):47–55.

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

The RCFPs (including DsRed-Express and DsRed-Express2) are covered by one or more of the following U.S. Patent Nos. 7,166,444; 7,157,565; 7,217,789; 7,338,784; 7,338,783; 7,537,915 6,969,597, 7,150,979 and 7,442,522.

Clontech has a license to sell products containing WPRE, under the terms described below. Any use of WPRE outside of Clontech's product or the product's intended use, requires a license as detailed below. Before using the product containing WPRE, please read the following license agreement. If you do not agree to be bound by its terms, contact Clontech within 10 days for authorization to return the unused product containing WPRE and to receive a full credit.

Patents: The WPRE technology is covered by one or more of the following U.S. Patents and corresponding patent claims outside the U.S.: 6,136,597; 6,284,469; 6,312,912; 6,287,814, issued to The Salk Institute for Biological Studies.

Individual License Agreement: Clontech grants you a non-exclusive license to use the enclosed product containing WPRE in its entirety for its intended use. The product is being transferred to you in furtherance of, and reliance on, such license. Any use of WPRE outside of Clontech's product or the product's intended use, requires a license from the Salk Institute for Biological Studies.

Termination of License: This license agreement is effective until terminated. You may terminate it at any time by destroying all products containing WPRE in your control. It will also terminate automatically if you fail to comply with the terms and conditions of the license agreement. You shall, upon termination of the license agreement, destroy all products containing WPRE in your control, and so notify Clontech in writing. This License shall be governed in its interpretation and enforcement by the laws of the State of California.

Contact for WPRE Licensing: The Salk Institute for Biological Studies, 10010 North Torrey Pines Road, La Jolla, CA 92037, Attn.: Office of Technology Management, Phone: 858.453.4100 ext. 1275, Fax: 858.546.8093.

This product and its use are the subject of U.S. Pat. No. 6,682,907. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot disclose information, sell or otherwise transfer this product, its components or materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for any commercial purposes. If the buyer is not willing to accept the limitations of this limited use statement, Clontech is willing to accept return of the product with a full refund. For information on purchasing a license to the DNA-Flap technology for purposes other than research, contact the Transfer of Technology Office, Institut Pasteur, 28 rue du Docteur Roux, 75 724 Paris Cedex 15 (www.pasteur.fr).

Living Colors Fluorescent Protein Products:

Not-For-Profit Entities: Orders may be placed in the normal manner by contacting your local representative or Clontech Customer Service at 650.919.7300. At its discretion, Clontech grants Not-For-Profit Entities a non-exclusive, personal, limited license to use this product for non-commercial life science research use only. Such license specifically excludes the right to sell or otherwise transfer this product, its components or derivatives thereof to third parties. No modifications to the protein coding sequence may be made without express written permission from Clontech. Any other use of this product requires a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com.

For-Profit Entities wishing to use this product are required to obtain a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com.

DH5 α TM and DH10BTM are trademarks of Life Technologies Corporation.

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc., unless noted otherwise. Clontech is a Takara Bio Company. ©2010 Clontech Laboratories, Inc.