

## Table of Contents

I. Introduction.....	1
II. Required Materials.....	1
III. Staining Protocol (35 mm Dish Format).....	1

## List of Tables

Table 1. Preparing X-Gal Staining Mix.....	2
--	---

## I. Introduction

The **Beta-Galactosidase Staining Kit** (Cat. No. 631780) provides a complete set of reagents for X-gal detection of beta-galactosidase expression in mammalian cells.

## II. Required Materials

### Supplied in Kit:

- X-Gal Solution (20 mg/ml)
- Staining Solution 1\*
- Staining Solution 2\*
- Staining Solution 3
- 10X Fixing Buffer\*
- 10X PBS

**\*CAUTION:** Due to their toxic properties, wear gloves and take appropriate precautions (see MSDS) when using Fixing Buffer (Section III, Steps 2–5) and Staining Solutions 1 and 2 (Section III, Steps 4–8).

### Not Supplied:

- Deionized H<sub>2</sub>O
- 37°C incubator
- Light microscope
- 35 mm cell culture dishes

## III. Staining Protocol (35 mm Dish Format)

1. Gently wash the cells with 2 ml of 1X PBS in a 35 mm\* cell culture dish.

\* Scale up as needed if you are using a larger dish.

**NOTE:** Washing too vigorously can cause detachment of less adherent cells.

2. Add 2 ml of 1X Fixing Buffer\* to the dish.

\* Prepare 1X Fixing Buffer by diluting 200 µl of 10X Fixing Buffer 1:10 with deionized H<sub>2</sub>O.

## Beta-Galactosidase Staining Kit Protocol-At-A-Glance

3. Incubate 5–10 minutes at room temperature.

**NOTE:** Longer fixation times can result in enzyme inactivation and loss of signal.

4. While the cells are fixing, prepare the appropriate amount of X-Gal Staining Mix (Table I).

**Table 1. Preparing X-Gal Staining Mix**

Component	1 Dish	6 Dishes
X-Gal Solution	100 µl	600 µl
Staining Solution 1	20 µl	120 µl
Staining Solution 2	20 µl	120 µl
Staining Solution 3	20 µl	120 µl
1X PBS	1.84 ml	11.04 ml
Total Volume	2 ml	12 ml

**NOTE:** If there is a visible precipitate after thawing the staining solutions, warm to 37°C and vortex until dissolved.

5. Wash the dish twice with 2 ml of 1X PBS.

**NOTE:** Thorough washing is necessary to remove the Fixing Buffer in order to prevent inhibition of the enzyme reaction.

6. Add 2 ml of the X-Gal Staining Mix to the dish.
7. Incubate the dish at 37°C for 1 hr to overnight.
8. Assay for development of blue color using light microscopy.

Contact Us For Assistance	
Customer Service/Ordering	Technical Support
Telephone: 800.662.2566 (toll-free)	Telephone: 800.662.2566 (toll-free)
Fax: 800.424.1350 (toll-free)	Fax: 650.424.1064
Web: <a href="http://www.clontech.com">www.clontech.com</a>	Web: <a href="http://www.clontech.com">www.clontech.com</a>
E-mail: <a href="mailto:orders@clontech.com">orders@clontech.com</a>	E-mail: <a href="mailto:tech@clontech.com">tech@clontech.com</a>

### Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without prior written approval of Clontech Laboratories, Inc.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product's web page at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

Clontech and the Clontech logo are trademarks of Clontech Laboratories, Inc. All other marks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech is a Takara Bio Company. ©2012 Clontech Laboratories, Inc.

This document has been reviewed and approved by the Clontech Quality Assurance Department.